MAS-100 CG Ex[®]

Air Monitoring for pressurised gases: Air / Nitrogen / Carbondioxyd / Oxygen / Argon

MAS-100 CG Ex®

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MAS-100 CG Ex®

The MAS-100 CG Ex® has been developed to control the microbiological contamination in compressed gases. The sampling method is based on the approved standard impaction principle, but under pressure. Standard 90-100mm Petri dishes can be used. The most widely used gases in the industry are pre-programmed. All instruments are delivered as explosion-proof versions.

Unique: The micro organisms are collected under the same pressure as in the gas-distribution system. After sampling, the automatic decompression cycle is performed.

New: Pre-programmed also for Oxygen (O₂)

Application field

Our product has been designed for compressed gases used in the pharmaceutical and food industries and other critical fields. Compressed gases may be connected directly to the instrument from 1 up to 10 bars. The instrument is easy to handle, works absolutely automatically and is operated by a rechargeable battery. It is possible to connect it to a PC.

Function al principle

The MAS-100 CG Ex® is an air sampler that is based on the impaction principle. The resulting airflow is directed onto a standard Petri dish containing agar. After the collection cycle, the Petri dish needs to be incubated and the colonies can be counted afterwards.

Advantages

Standard explosion proof version
Integrated pressure and volume sensors
Uses standard 90-100mm Petri dishes
Dialogue software
PC interface

Technical specifications

Airflow 100 or 50 litres/minute at 1 –10 bar
Rechargeable battery
Autonomy approx. 3.5h
Explosion-proof, powder-coated aluminium case
Sampling head sterilisable and with pressure certificate
Pre-programmed gases: Air, N₂, CO₂, Ar, O₂
ATEX certified

Dimensions

without sampling head: 33 x 23 x 11 cm
with sampling head: 49 x 32 x 11 cm
Weight with sampling head: Approx. 12.5 kg
Dimensions of transport trolley: L 62 x H 42 x W 26 cm

Firmware

5 freely programmable volumes in the range of 1-2000 litres
Interactive menu
Real-time clock with date and calibration reminder
PC interface

Accessories

Transport trolley (standard)
Additional sampling head (optional)