

ADAPTER FOR GLASS PETRI DISHES FOR VIABLE AIR SAMPLERS MAS-100 /NT/EX AND MAS-100 ECO

PRODUCT INFORMATION

The adapter is padded with rubber discs to hold glass Petri dishes in place without damaging them. It permits holding glass Petri dishes in place on these viable air samplers: MAS-100, MAS-100 NT, MAS-100 NT Ex and MAS-100 Eco.


The geometry of the glass Petri dish has to be within these ranges:

Outer diameter: 90 mm - 93 mm

Max. height: 18 mm (without lid)

INSTALLATION INSTRUCTION

1. Remove perforated lid from the sampling head.
2. Using the Allen key, unscrew the three screws holding the blue standard jaws (below right). Take care not to lose the springs. The screws and springs are used to fix the special jaws (below left) included in this set.
3. Replace the standard jaws with the special jaws included in the set. The rubber pads point to the center of the plate holder.
4. Adjust the three jaws with the Allen key such that the Petri dish is snugly held in place. Do not tighten the jaws too much.
5. Adjust the three jaws such that the Petri dish is centered in the plate holder.

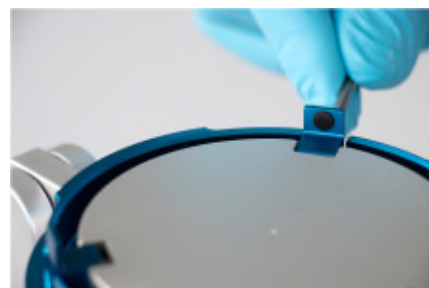

ORDER INFORMATION

Name	Merck Article number	MBV Article Number
Adapter for glas Petri dish	1.09082.0001	06.6098.01
Clamp blue for standard Petri dish with screws and springs	1193830001	06.6099.01